FLEXIBLE-ARM[™] LIGHT INSTRUCTIONS

SPECIFICATIONS:

Electrical:	150 VAC, 60Hz, UL Listed
Light:	35 watt Halogen bulb
Plug:	Polarized
Arm length:	Cat. No. 3095 - 21"
U U	Cat No. 3096 - 31"
Weight:	8 lbs.

FEATURES

- Base 7" diameter, 7lb.
- 35 watt bright Halogen white light
- Clamping unit for vertical or horizontal Positioning
- Switch in cord
 Flexible-Arm [™] articulates at 3 points
- · Fingertip bar for raising, lowering and
- pivoting with ease
- Spring counterbalance for stability
- Vented shade with rolled edges
- · Safety-glass with UV filter

- BASE ASSEMBLY (see Figure 1) 1. Remove "Screw I" from base "Stem II".
- 2. Insert the lamp frame into base "Stem II",
- align screw holes.
- 3. Screw in "Screw I" and finger tighten.

CLAMPING UNIT ASSEMBLY

- 1. Insert the lamp frame into hole in top of clamp assembly.
- 2. Loosen clamp with lever on bottom of assembly.
- 3. Position clamping mechanism and tighten firmly using lever on bottom of assembly.

BULB REPLACEMENT (see Figure 2) 1. Turn off lamp and allow shade and

- accessories to cool.
- 2. Turn the lampshade upward.
- 3. Remove CENTER screw (using a Phillips screwdriver) and clip from protective glass cover.
- 4. Pull out failed bulb and replace with
- halogen bulb. (Control Company Cat. No. 3098) 5. Replace protective glass and clip and screw in place.

NOTE: Do not remove screws holding fingertip bar in place.

WARRANTY, SERVICE, OR RECALIBRATION

For warranty, service, or recalibration, contact:


CONTROL COMPANY

4455 Rex Road Friendswood. Texas 77546 USA Ph. 281-482-1714 Fax 281-482-9448 E-mail sales@control3.com www.control3.com

Control Company is ISO 9001 Quality-Certified by DNV and ISO 17025 accredited as a Calibration Laboratory by A2LA.

> Cat. No. 3095/3096 Flexible-Arm™ Light Instructions 92-3095-10 Rev. 1 071504

Traceable® is a registered trademark of Control Company

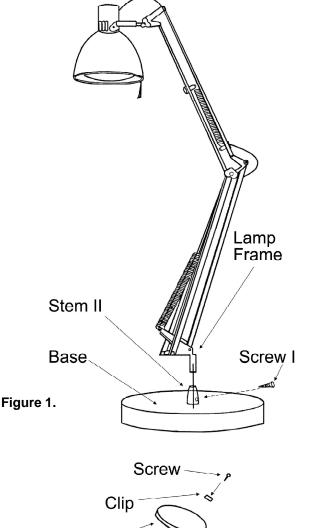

Protective Glass - Halogen Bulb Shade Assembly Shade -

Figure 2.

© 2004 🕑 Control Company