

SPECIFICATIONS


Display: 1/4"-high, 4-digit
Time Range: 19 hrs to 59 mins
Resolution: 1 minute
Accuracy: 0.01%
Size: 2 x 1 1/8 x 3/8 inches
Weight: 1/2 ounce
Features: Key-chain, Count down/Countup Timing

Controls:

H - Hour button.

M - Minute button.

S/S - Start/Stop button.

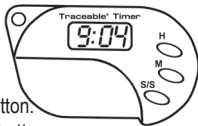

OPERATION

Clearing The Display:

To clear the display to 0:00 press the H and M buttons simultaneously. The unit will beep 2 (two) times when the display is cleared.

COUNTDOWN TIMER/ALARM OPERATION

1. Clear the display to 0:00.
2. Press H button to set number of hours desired. Setting will advance 1 hour each time button is pressed. Press and hold to advance rapidly.
3. Press M to set number of minutes

desired. Setting will advance 1 minute each time button is pressed. Press and hold to advance rapidly.

4. Once desired time is displayed, press the S/S button to start the countdown operation.
5. At zero the alarm will sound for 1 minute and stop automatically.
6. Pressing the H and M buttons simultaneously silences the alarm and clears the display to zero.
7. Press the S/S button to manually silence the alarm and to also begin countdown timing again from the time set.

Note: The blinking ":" symbol in the display indicates timer is running.

Count-up/Stopwatch Operation

1. Clear the display to 0:00.
2. Press the S/S button to start count-up/stopwatch operation.
3. Press S/S button to stop count-up/stopwatch operation.

Note: When timing capacity reaches 19 min. and 59 sec. the timer automatically rolls over to zero and continues timing.

TIME OUT

If a pause in countdown or count-up operation is required press the

SS button to “freeze” display at current reading (time-out). To restart operation, press S/S button again. An unlimited number of time-outs may be taken.

OPERATIONAL DIFFICULTIES

If this unit does not function properly for any reason, replace the battery with a new high-quality battery (see “Battery Replacement” section). Low battery power can occasionally cause any number of “apparent” operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

BATTERY REPLACEMENT

Erratic readings, a faint display, no display, or a flashing display are all indicators that the battery must be replaced. Slide battery cover off the backside of unit in direction indicated by arrow. Replace exhausted battery with a new 1.5-volt Battery with the positive side up. Replace battery cover. Replacement battery Cat. No. 1104.

WARRANTY, SERVICE, OR RECALIBRATION

For warranty, service, or recalibration, contact:

TRACEABLE® PRODUCTS

12554 Old Galveston Rd. Suite B230

Webster, Texas 77598 USA

Ph. 281 482-1714 • Fax 281 482-9448

E-mail support@traceable.com

www.traceable.com

Traceable® Products is ISO 9001:2015 Quality-Certified by DNV and ISO/IEC 17025:2005 accredited as a Calibration Laboratory by A2LA.

Cat. No. 1060

Traceable® is a registered trademark of Cole-Parmer.

©2018 Traceable® Products.
92-1060-00 Rev. 3 061418

TRACEABLE® KEY-CHAIN TIMER INSTRUCTIONS