


# Download TraceableLIVE mobile app via Apple App Store or Google Play

1

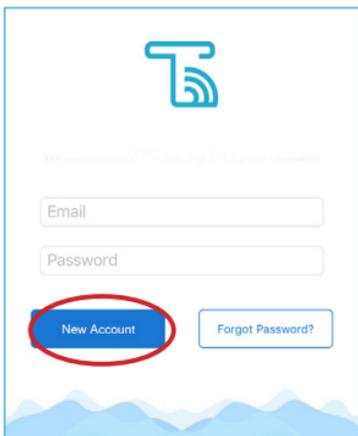


Register Account

# Add Device to Account

2

Open the TraceableLIVE app:

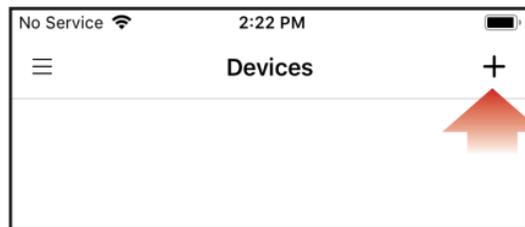

Tap the + icon to add a device.

- Permission must be given to TraceableLIVE app to access your mobile device's camera.
- Using your mobile device's camera hover over QR code on back of unit (see right).
- App will automatically fill in device information.



# WiFi Setup

3



Press/HOLD WiFi button "AP" will appear on display.

Press WiFi button again, "AP URIF" will appear on display.

When "AP rEAdY" appears, proceed to next step.


# Use Internet enabled device to connect to WIFI

4



Select Network: CC6500-0013 see example network

Example Network: CC6500-0013

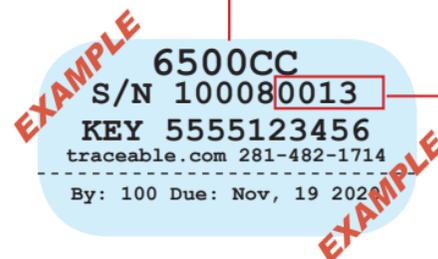

Open Browser, type  
192.168.1.1 in address bar

5

Select local network and enter password.

Security type WPA2 recommended or  
contact your network admin.



**CONGRATULATIONS!**  
Your device will restart.

**Still having trouble?**  
Please give us a call  
281-482-1714 or email  
[support@traceable.com](mailto:support@traceable.com)

**TraceableLIVE®**  
**QUICK START**  
**GUIDE**