


Go to traceablelive.com to create a new account.

1

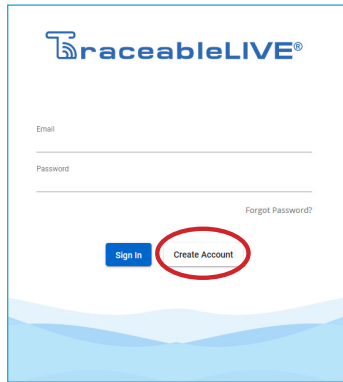

TraceableLIVE®

Email

Password

[Forgot Password?](#)

[Sign In](#) [Create Account](#)


Register Account

THEN Download TraceableLIVE mobile app via Apple App Store or Google Play.

Add Device to Account

2

Open TraceableLIVE app on mobile or desktop and Log in to account:

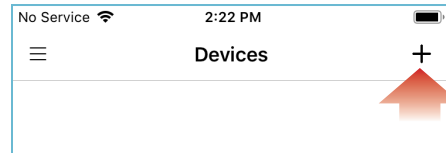

Tap the **+** icon to add a device.

- Permission must be given to TraceableLIVE app to access your mobile device's camera.
- Using your mobile device's camera hover over QR code on back of unit (see right).
- App will automatically fill in device information.




9210382

Note: QR code only works with a mobile device.

Ethernet Cable

3

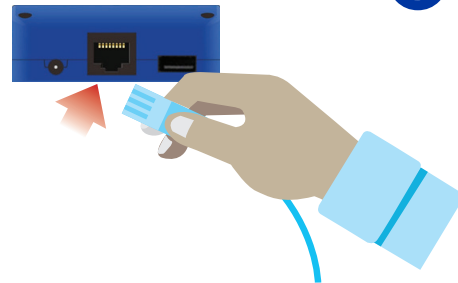

INSERT ETHERNET CABLE

- On back of unit, locate Ethernet cable jack.
- Line up the cable connector with the jack and push into jack until plug clicks into place.

INSERT AC ADAPTOR

- AC adaptor must be plugged in as the primary power source.

INSERT PROBE

- Insert probe into USB jack.

CONGRATULATIONS!

4

Device will start to post data if AC adaptor and Ethernet cable are plugged in, and an active link to server is established.

TraceableLIVE® QUICK START GUIDE



DEVICE ONLY

**Still having trouble?
Please give us a call
281-482-1714 or email
support@traceable.com**