TURNING THE UNIT ON/OFF To turn the unit on, press the ON button.	INCHES OR MILLIMETERS Once the unit has been turned on, each press of the INCH/MM button will toggle the display	ZERO ADJUSTING Clean the measuring surfaces by rubbing a clean piece of paper between the caliper surfaces	NOTE The graduated scale on the caliper body is for approxim reference only, the exact
To turn the unit off, press the OFF button.	between inches and millimeters.	several times.	measurement, accurate to ±0.2mm/0.01", is shown on
	While taking a measurement, press the INCH/MM button at any time to covert from one unit of	Push the measuring surfaces together.	LCD display.
	measure to the other.	Reset the display to zero by pressing the 0 button.	

ALL OPERATIONAL DIFFICULTIES

If this unit does not function properly for any reason, replace the battery with a new high-quality battery (see "Battery Replacement" section). Low battery power can occasionally cause any number of "apparent" operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

BATTERY REPLACEMENT

Erratic readings, a faint display, or no display are all indications that the battery must be replaced. Slide the battery cover off (away from the display). Remove the exhausted battery and replace with a new 1.5 volt silver-oxide battery. Equivalent battery replacements are: RAYOVAC RW42, DURACELL D357, GPA76, and EVEREADY 357. Make certain to insert the new battery with the positive (+) side up. Replace the battery cover.

WARRANTY, SERVICE, OR RECALIBRATION

For warranty, service, or recalibration contact:

TRACEABLE® PRODUCTS 12554 Old Galveston Rd. Suite B230

Webster, Texas 77598 USA
Ph. 281 482-1714 • Fax 281 482-9448
E-mail support@traceable.com • www.traceable.com

Traceable® Products are ISO 9001:2018 Quality-Certified by DNV and ISO/IEC 17025:2017 accredited as a Calibration Laboratory by A2LA.

Traceable® is a registered trademark of Cole-Parmer.

©2020 Traceable® Products, 92-3417-10 Rev. 5 042220

TRACEABLE® CARBON-FIBER CALIPER INSTRUCTIONS