

SPECIFICATIONS

File length: 4 inches

File width: ¼" half round

Handle length: 4 inches

Length of entire unit: 8 inches

Glassware Rescuer™ is constructed using real diamonds; the hardest material known. Use it to repair all beakers, flasks, and lab glassware, as well as most hard materials.

OPERATIONS

1. Moisten Glassware Rescuer™ with water.
2. With round side down, rub back and forth in the recessed edges of the broken object.
3. With flat side down, rub back and forth on the sharp raised corners of the broken object.

NOTE: Only a few strokes should be needed to smooth most surfaces.

CAUTION: Eye protection and gloves should be worn. Caution should be used when repairing all glassware.

**WARRANTY, SERVICE, OR
RECALIBRATION**

For warranty, service, or recalibration, contact:

TRACEABLE® PRODUCTS

12554 Old Galveston Rd. Suite B230

Webster, Texas 77598 USA

Ph. 281 482-1714 • Fax 281 482-9448

E-mail support@traceable.com

www.traceable.com

Traceable® Products are ISO 9001:2018 Quality-
Certified by DNV and ISO/IEC 17025:2017
accredited as a Calibration Laboratory by A2LA.

©2020 Traceable® Products. 92-3315-00 Rev. 4 042420

**GLASSWARE
RESCUER™
INSTRUCTIONS**