
SPECIFICATIONS


OPERATION

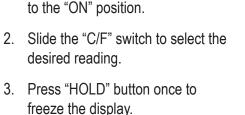

4. Press the "HOLD" button again to

resume measurement.

Slide the power "ON/OFF" switch

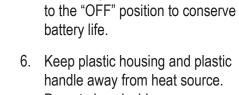


battery life. Keep plastic housing and plastic handle away from heat source. Do not place inside oven or expose to open flame.

Do not immerse probe handle or

cable in liquid.

Slide the power "ON/OFF" switch

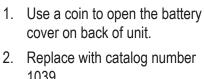

cover on back of unit. 2. Replace with catalog number 1039.

BATTERY REPLACEMENT

WARRANTY, SERVICE, OR RECALIBRATION

For warranty, service, or recalibration, contact:

TRACEABLE® PRODUCTS

12554 Old Galveston Rd. Suite B230 Webster, Texas 77598 USA Ph. 281 482-1714 • Fax 281 482-9448 E-mail support@traceable.com

Traceable® Products are ISO 9001:2018 Quality-Certified by DNV and ISO/IEC 17025:2017 accredited as a Calibration Laboratory by A2LA. TRACEABLE®
MINITHERMOMETER
INSTRUCTIONS

Traceable® is a registered trademark of Cole-Parmer.

©2020 Traceable® Products. 92-4146-00 Rev. 3 050120