

SPECIFICATIONS

Range:

–58 to 302°F (–50 to 150°C)

Resolution

0.1° from –20.0 to 100°C

Accuracy:

±1°C from –20 to 200°C

OPERATION

1. Slide the power “ON/OFF” switch to the “ON” position.
2. Slide the “C/F” switch to select the desired reading.
3. Press “HOLD” button once to freeze the display.
4. Press the “HOLD” button again to resume measurement.
5. Slide the power “ON/OFF” switch to the “OFF” position to conserve battery life.
6. Keep plastic housing and plastic handle away from heat source. Do not place inside oven or expose to open flame.
7. Do not immerse probe handle or cable in liquid.

BATTERY REPLACEMENT

1. Use a coin to open the battery cover on back of unit.
2. Replace with catalog number 1039.

**WARRANTY, SERVICE, OR
RECALIBRATION**

For warranty, service, or recalibration, contact:

TRACEABLE® PRODUCTS

12554 Old Galveston Rd. Suite B230

Webster, Texas 77598 USA

Ph. 281 482-1714 • Fax 281 482-9448

E-mail support@traceable.com

www.traceable.com

Traceable® Products are ISO 9001:2015 Quality-
Certified by DNV and ISO/IEC 17025:2017
accredited as a Calibration Laboratory by A2LA.

Traceable® is a registered trademark of Cole-Parmer.

©2020 Traceable® Products. 92-4146-00 Rev. 4 072524

**TRACEABLE®
MINI-
THERMOMETER
INSTRUCTIONS**