

SPECIFICATIONS

Probe Range: -58.0 to 158.0°F / -50.0 to 70.0°C

Resolution: 0.1°

Update Rate: 10 seconds

Battery: AAA (1.5V)

INSTALLATION

1. Use sensor holder to fix external sensor at desired external space.
2. Plug external sensor into unit.
3. Pull out the battery insulation strip.
4. Unit will perform a self-test and display readings normally after 3 seconds.
5. Place unit away from direct sunlight, rain or extreme heat.
6. Peel off display protective sheet.

DISPLAYING °F OR °C

To display temperature readings in Fahrenheit or Celsius, press °C/°F button located on front of unit.

MINIMUM AND MAXIMUM MEMORY

1. Press Max/Min to display maximum measured value.  will appear on display. Press button once more to display minimum measured value.  icon would appear. Press Max/Min once more to return to current temperature.

2. Press and hold [Max/Min] until two bars -- are displayed then release to reset memory.

TEMPERATURE ALARM CONFIGURATION

To enter configure alarm mode press and hold, [ALERT/SET] () will begin to flash on display.

1. Press °C/°F button to set HIGH alarm setting.
2. Press [ALERT/SET] setting will flash.
3. Press °C/°F button to set LOW alarm setting.
4. Press [ALERT/SET] to finish configuration.

Notes: When temperate is OUTSIDE configured limit an audible alarm will beep for 5 seconds.

Press and hold Hold [°C/°F] to increase the value automatically.

TO TURN ON/OFF ALARMS

1. To deactivate the alarm, press [ALERT/SET] to put the screen in off position.
2. To switch on the alarm: Press [ALERT/SET] to put the screen in on position.
3. The screen ALE OFF or ALE ON will return to the normal screen automatically.

WARRANTY, SERVICE, OR RECALIBRATION

For warranty, service, or recalibration, contact:

TRACEABLE® PRODUCTS

12554 Old Galveston Rd. Suite B230

Webster, Texas 77598 USA

Ph. 281 482-1714 • Fax 281 482-9448

E-mail support@traceable.com

www.traceable.com

Traceable® Products are ISO 9001:2015 Quality-
Certified by DNV and ISO/IEC 17025:2017
accredited as a Calibration Laboratory by A2LA.

Cat. No. 4126

©2020 92-4126-00 Rev. 8 100325

**TRACEABLE®
BIG-DIGIT
MEMORY
THERMOMETER
INSTRUCTIONS**